

Technical Datasheet

EZAssay[™] Reactive oxygen species (ROS) Assay Kit

Product Code: CCK078

1. Introduction:

Reactive oxygen species (ROS) are produced inside the cell as a metabolic by product of oxidation process. When ROS production goes above threshold level, it causes oxidative stress and damages the cellular integrity. ROS contribute to various cell signaling process and induces various disease conditions like Parkinson's and Alzheimer's disease, aging, cancer etc. Presence of ROS in cellular environment can be detected using cell permeable fluorogenic dye. Dye is sensitive towards hydroxyl, peroxyl and another ROS activity within the cell. Fluorogenic dye can be detected by using a cytation, spectrofluorimeter or flow cytometer. The excitation spectra is 485nm and while emission spectra is 528nm.

2. About the Assay:

EZAssayTM ROS assay Kit is designed for determination of ROS activity within a cell using a cell permeable fluorogenic dye. Once fluorogenic dye diffuses inside the cells, esterases deacetylate it to a non-fluorescent compound. In presence of ROS, the non-fluorescent compound immediately gets oxidized to a highly fluorescent product, which is easily detectable at excitation wavelength of 485nm and Emission wavelength of 528nm. The fluorescent intensity produced is directly proportional to ROS level in the cells.

3. Applications:

• Cell metabolism: Comparative quantification of changes in cellular activity of with or without treatment of any compound or drug.

• Cell cytotoxicity: Evaluation of effect of inhibitors or inducers of apoptosis.

• Drug discovery: High-throughput screening of various anticancer, anti-inflammatory drugs.

4. Kit contents:

Reagents supplied in the kit are sufficient for 200 assays (96microwell plates-2 Nos.).

Code	Description	Quantity	Storage
CCK078(A)	ROS staining reagent	2 x 20 µ1	-30°C to -10°C
CCK078(B)	Active oxygen positive control (Tert-butyl hydrogen peroxide)	0.5 ml	15°C to 30°C
CCK078(C)	Washing buffer	100.0 ml	15°C to 30°C

5. Materials required but not provided in the kit:

- Cells in appropriate medium without phenol red.
- Adjustable pipettes and a repeat pipettor.
- Transparent flat-bottom black 96-well microtiter plate for culturing the cells.
- 96-well plate reader capable of measuring the fluorescence at 450nm and >600nm.

6. General guidelines:

It is important to optimize experimental factors like cell density, incubation time, media composition and concentration of the agents under investigation prior to use of ROS Assay Kit.

6.1 Assay controls

Include appropriate assay controls i.e.

- 1. Cell control (medium with cells but without the experimental drug /compound and fluorogenic dye).
- 2. Vehicle control (medium containing the experimental drug or compound with cells).
- Positive control (medium with cells + 5 to 20µl of CCK078(B) per well)
- 4. Negative control Medium (without cells & experimental drugs) and fluorogenic dye.

6.2 Accuracy

• Perform the assay in triplicates or more to obtain statistically significant data.

• Accuracy of the assay depends on pipetting skills of the personnel. Inappropriate addition and mixing practices may result in erroneous and false-positive or false-negative results.

• Use of a repeating pipettor is recommended to pipette reagents. This saves time and helps to maintain more precise incubation times.

Pipette tip should be equilibrated with the reagent before use. This is carried out by slowly filling up the tip with reagent and gently expelling the contents several times.
Care should be taken so that no bubbles are introduced into the wells during pipetting or mixing of the reagents.

6.3 Culture Medium

Phenol red may interfere with fluorogenic dye. Therefore, the cell culture media used for this assay should not contain phenol red.

6.4 Temperature

Temperature affects the performance of the assay because of its effect on enzymatic reaction rates. It is crucial to run the assay at a uniform temperature to ensure reproducibility across a single plate or among stacks of several plates. Since fluorescence readings are measured at room temperature, it is important to ensure adequate equilibration of assay plates after removal from a 37°C incubator to avoid differential temperature gradients. Stacking large numbers of assay plates in close proximity should be avoided to ensure complete temperature equilibration.

6.5 Measurement of absorbance

Fluorescence can be read at an excitation wavelength of 485nm and emission at 528nm.

7. Directions for use

Users are advised to review entire procedure before starting the assay

7.1Preparation of working CCK078(C)1ml of CCK078(C) + 9ml of cell culture grade water.

7.2 Preparation of cells

Always use freshly harvested cells for assay. Seed the cells in a cell culture flask or dish in an amount appropriate for the assay and incubate at 37°C in a 5% CO2 environment. Allow the cells to grow for up to 24 hours or till confluency reaches. Harvest the cells and use for assay. *Note: The quantity of the cell suspension to be seeded in the medium depends upon doubling time of individual cell lines and seeding density to be used in the assay.*

7.3 Assay procedure

- 1. Harvest the cells as explained in section 7.2.
- 2. Adjust appropriate cell density as per the user requirement and type of cells being used
- 3. Seed 100µl cells/well in a 96 well black well plate with transparent bottom.
- 4. After 24 hours of incubation, discard the spent medium, wash once with 1X CCK078(C).
- 5. Add 100µl of 1X CCK078(C) or desired medium.
- Add test molecule or a positive control i.e. CCK078(B) in the test wells, (concentration of test molecule should be as per user's requirement) and incubate the plate at 37°C with 5% CO2.
- For positive control, add 5 to 20 μl of CCK078(B) and incubate the plate for 30 minutes at 37°C with 5% CO₂.
- Thaw one bottle of CCK078(A) and equilibrate at room temperature just before use.
 Note: Reagent in one bottle is sufficient for one 96-well plate.
- Take 12.5 μl of CCK078(A) and add in 10ml of medium (without serum) and make a homogenous solution.

Note: CCK078(A) is light sensitive hence, make sure that experiments are carried out in dark.

- 10. Remove the test molecule/positive control from wells
- 11. Wash the cells once with working solution of CCK078(C).
- Add 100µl of homogenous solution of fluorogenic dye to each well except in cell control and mix the dye uniformly by rotating the plate (clockwise and anti-clockwise).
- 13. Wrap the plate with aluminum foil to avoid exposure to light.
- 14. Incubate the plate at 37°C for 20 to 30 minutes.
- 15. After incubation, measure the fluorescence at Ex/Em = 485 nm/528 nm.
- Subtract the blank (cell control) value to determine the reading due to fluorescence. Calculate the mean.
- 17. Plot a graph of normalized mean fluorescence versus cell density.



Fig. 1: CHO cells were seeded at different densities per well in a black well clear bottom 96 well plate. Each cell density was treated with 5 μ L of CCK078(B) for 30 min at 37°C, 5% CO2. After 30 min, the cells were washed once with 1X CCK078(C). ROS staining was performed and fluorescence intensity was measured using Cytation reader at excitation 485nm and emission 528nm.

8. Storage and shelf life

Store CCK078(A) dye at -30°C to -10°C once received. Repeated freezing and thawing of CCK078(A) may result in loss in activity of the reagent and increased background absorbance.
Use before expiry date.

9. Advantages

Time saving: Fluorescence can be measured without involving solubilization with organic solvent. **Easy reagent preparation:** Ready to mix reagents offer ease of reagent preparation.

Reproducibility: Entire assay can be performed in a single plate. Cells and reagents need not be transferred. This facilitates reproducibility of the results. **Sensitivity and accuracy:** ROS staining reagent strongly correlate with the metabolic activity of the cells. This allows use of low cell densities.

Safety: No radioisotopes are involved. Fast: Use of black transparent flat bottom 96 well plates allow the processing of large number of samples. Flexibility: ROS staining reagent works on adherent as well as suspension cell lines.

10. Troubleshooting points

Use the following troubleshooting guidelines for technical assistance

Problem	Cause	Solution
Color change in ROS staining reagent	Exposure of the reagent to light.	Wrap the reagent bottles and culture plates with aluminum foil.
Inappropriate readings	Improper selection of the filter.	Choose appropriate filters.
	Cross contamination due to inaccurate pipetting technique or inaccurate equipment.	Perform the assay using automated electronic pipettes for seeding the cell suspension and adding the reagents.
Very high fluorescence values	Microbial contamination	Discard. Repeat the assay with new media and reagents.
	Too much signal in ROS staining reagent signal due to high cell densities.	Repeat the assay with reduced cell densities.
Random fluorescence values/ poor consistency of replicates	Test compound under study is responsible for improper response of the cells to ROS staining reagent.	Refer to the pharmacological properties of the compound.
	Microbial contamination	Discard. Repeat the assay with new media and reagents.

Disclaimer:

Revision : 06/2025

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia[™] publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia[™] Laboratories Pvt. Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.



Plot No. C40, Road No. 21Y, MIDC, Wagle Industrial Area, Thane (West) 400604, Maharashtra, India.Tel No.022-69034800 Email: atc@himedialabs.com Website: www.himedialabs.com.