

## HiIndicator™ pH papers

LA310, LA312, LA315, LA318, LA321, LA323, LA334, LA335.

The convenience of using HiIndicator papers for the rapid determination of pH values has led to many applications in laboratories and industry. These pH papers are made with special indicator dyes that change color at specified pH value.

Somewhat uneven colour of the strips is of no consequence. The colour obtained on use is indicative of the correct pH.

**Application :** Analytical chemistry, biology & various laboratories and industries etc.

Product Name	Product Code	Description	pH Range
HiIndicator™ pH papers.	LA310	HiIndicator pH paper	2.00 – 10.50
	LA312	HiIndicator pH paper	3.50 – 6.00
	LA315	HiIndicator pH paper	3.80 – 5.30
	LA318	HiIndicator pH paper	5.00 – 7.50
	LA321	HiIndicator pH paper	6.50 – 9.00
	LA323	HiIndicator pH paper	8.00 – 10.50
	LA334	HiIndicator pH paper	2.00 - 4.50
	LA335	HiIndicator pH paper	1.00 - 14.00

**Direction for use :** Tear off strip of indicator paper and insert it for a few seconds into the solution to be tested. With highly viscous or stained liquids and with suspensions, drip the substance onto the indicator paper. Compare the wet paper with the colour scale. For papers where liquids are dripped, compare the reverse side. Possible discolouration of the dry new papers may be caused by their high sensitivity. This does not impair the efficacy of the Indicator papers for pH determinations.

The so-called indicator error may occur with very weakly buffered or unbuffered solution and can be compensated for up to a point in the following manner. : - The strip can be made to adhere to the inner wall of the a test tube, which can then filled to the upper edge of the paper with the fluid to be tested. After 1/2 to 1 minute, the colour of the paper may be compared with the scale through the glass of test tube.

**Product Features:**

- Instant pH readings.
- Accurate for a wide range of routine pH testing.
- Convenient and portable for field use.
- Pack Size : 1 pack-200 Nos.

**Disclaimer :**

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.

---

HiMedia Laboratories Pvt. Ltd. A-516, Swastik Disha Business Park, Via Vadhani Ind. Est., LBS Marg, Mumbai-400086, India.  
Customer care No.: 022-6147 1919 Email: [techhelp@himedialabs.com](mailto:techhelp@himedialabs.com) Website: [www.himedialabs.com](http://www.himedialabs.com)