

MB522

HiPurA® Sperm Genomic DNA Purification Kit

Kit Contents

Product Code	Reagents provided	MB522		
		20 Preps	50 Preps	250 Preps
ML116	Resuspension Solution (1X PBS)	4 ml	10 ml	50 ml
DS0038	Semen Wash Buffer Concentrate (SEW)	12 ml	30 ml	150 ml
DS0039	Sperm Lysis Buffer (SL)	4 ml	10 ml	50 ml
DS0010	Lysis Solution (C1)	6 ml	15 ml	75 ml
DS0012	Wash Solution Concentrate (WS)	10 ml	25 ml	125 ml
DS0040	Elution Buffer (ET) [10 mM Tris-Cl, pH 8.5]	6 ml	15 ml	75 ml
MB086	Proteinase K	10 mg	25 mg	125 mg
DS0003	RNase A Solution (20 mg/ml)	0.6 ml	1.5 ml	7.5 ml
DBCA03	HiElute Miniprep Spin Column (Capped) [in DBCA016 Collection Tube]	20 nos	50 nos	250 nos
DBCA016	Collection Tube (Uncapped), Polypropylene (2.0 ml)	20 nos	50 nos	250 nos
PW1139	Collection Tube, Polypropylene (2.0 ml)	40 nos	100 nos	2 X 250 nos

Intended Use

Recommended for isolation of DNA from sperm samples of human and animal origin.

Introduction

HiPurA® Sperm Genomic DNA Purification Kit provides a fast and easy method for purification of total DNA for reliable applications in PCR and Southern blotting technique. The DNA purification procedure using the miniprep spin column comprises of three steps viz. adsorption of DNA to the membrane, removal of residual contaminants and elution of pure genomic DNA. HiMedia's HiElute Miniprep Spin Column (Capped) format allows rapid processing of multiple samples. The columns have a high binding capacity and high quality genomic DNA is obtained from various species. The DNA obtained is compatible with downstream applications such as restriction endonuclease digestion, PCR and Southern blotting.

HiPurA® Sperm Genomic DNA Purification Kit

This kit simplifies isolation of DNA from sperm samples with spin column procedure. Following lysis, DNA binds to the silica gel membrane of the HiElute Miniprep Spin Column (Capped) to yield high purity DNA. Two rapid wash steps remove trace salts and protein contaminants resulting in the elution of high quality DNA in the Elution Buffer (ET) provided with the kit.

HiElute Miniprep Spin Column (Capped) [DBCA03]

HiElute Miniprep Spin Column (Capped) is based on the advanced silica binding principle presented in a microspin format. The system efficiently couples the reversible nucleic acid-binding properties of the advanced silica gel membrane and the speed plus versatility of spin column technology to yield high quantity of DNA. The use of spin column facilitates the binding, washing and elution steps thus enabling multiple samples to be processed simultaneously. This column eliminates the need for alcohol precipitation, expensive resins, and harmful organic compounds such as phenol and chloroform, otherwise employed in traditional DNA isolation techniques. DNA binds specifically to the advanced silica-gel membrane while contaminants pass through. PCR inhibitors such as divalent cations and proteins are completely removed in two efficient wash steps, leaving pure nucleic acid to be eluted in the buffer provided with the kit.

Elution

Elution with 200 µl of Elution Buffer (ET) or Molecular Biology Grade Water (ML024) will provide sufficient DNA to carry out multiple amplification reaction. Elution with volume less than 200 µl will increase the final DNA concentration, but will reduce the overall DNA yield.

Concentration, yield and purity of DNA

Spectrophotometric analysis and agarose gel electrophoresis will reveal the concentration and the purity of genomic DNA. Use Elution Buffer to dilute samples and to calibrate the spectrophotometer, measure the absorbance at 260 nm, 280 nm, and 320 nm using a quartz microcuvette. Absorbance readings at 260 nm should fall between 0.1 and 1.0. The 320 nm absorbance is used to correct for background absorbance. An absorbance of 1.0 at 260 nm corresponds to approximately 50 µg/ml of DNA. The $A_{260}-A_{320}/A_{280}-A_{320}$ ratio should be 1.6 –1.9. Purity is determined by calculating the ratio of absorbance at 260 nm to absorbance at 280 nm. DNA purified by HiPurA® Sperm Genomic DNA Purification Kit is free of protein and other contaminants that can inhibit PCR or other enzymatic reactions.

Concentration of DNA sample (µg/ml) = 50 x A_{260} x dilution factor.

Materials needed but not provided

- Incubator at 55°C
- 70°C water bath or heating block
- Tabletop Microcentrifuge (with rotor for 2.0 ml tubes)
- Ethanol (96 – 100%)
- Corex® (or any other suitable material) centrifuge tubes
- Lymphocyte Separation Medium (LSM) (Product Code: LS001) (Optional)
- Clean and dry bottle of appropriate capacity (for dilution of SEW Buffer) (**Refer point 5 of General Preparation Instructions**)
- Molecular Biology Grade Water (Product code: ML024)
- DTT (DL-Dithiothreitol) (MB070)

Storage

Store the HiPurA® Sperm Genomic DNA Purification Kit between 15-25°C except certain components as specified on each labels. Under recommended condition kit is stable for 1 year.

General Preparation Instructions

1. **Set the incubator at 55°C**
(For incubation of semen sample / semen pellet with Buffer SL to form semen lysate)
2. Preheat a water bath or heating block to 70°C.

3. **Thoroughly mix reagents**

Examine the reagents for precipitation. If any kit reagent forms a precipitate (other than enzymes), warm at 55-65°C until the precipitate dissolves and allow cooling to room temperature (15-25°C) before use.

4. **Reconstitute DTT (DL-Dithiothreitol) (not provided)**

Resuspend the DTT (DL-Dithiothreitol) powder in Molecular Biology Grade Water (Product code: ML024) to obtain a 1M stock solution.

Number of Preps	DTT (DL-Dithiothreitol)	Molecular Biology Grade Water
20	49 mg	0.32 ml
50	124 mg	0.80 ml
250	618 mg	4 ml

The product as supplied is stable at room temperature (15-25°C), upon reconstitution store at -20°C in dark (wrapped in an aluminium foil), as mentioned in storage instructions.

NOTE: The 1M DTT solution must be added directly to each sample preparation every time.

5. **Reconstitute Proteinase K (MB086)**

The HiPurA® Sperm Genomic DNA Purification Kit contains Proteinase K. Intensive research has shown that it is the optimal enzyme for use with the Lysis Solution provided in the kit. It is completely free of DNase and RNase activity. Proteinase K is the enzyme of choice for use with an SDS containing Lysis Solution. The specific activity of the Proteinase K is 33.5 units/mg dry weight.

Resuspend the Proteinase K (MB086) powder in Molecular Biology Grade Water (ML024) to obtain a 20 mg/ml stock solution.

Number of Preps	Proteinase K	Molecular Biology Grade Water
20	10 mg	0.5 ml
50	25 mg	1.25 ml
250	125 mg	6.25 ml

The product as supplied is stable at room temperature (15-25°C), upon reconstitution store at -20°C as mentioned in storage instructions.

NOTE: The Proteinase K solution must be added directly to each sample preparation every time. Do not combine the Proteinase K and Lysis Solution for storage.

6. **Preparation of Sperm Lysis Buffer (SL) (DS0039)**

To 100 µl of Sperm Lysis Buffer (SL), add 20 µl of the Proteinase K solution (20 mg/ml) (**Refer to step 5 of General Preparation Instructions**) and 8 µl of 1M DTT, just before use. Mix thoroughly by vortexing for 10-15 seconds.

NOTE: Sperm Lysis Buffer (SL) can be used effectively for semen as well as sperm sample.

7. **Dilute Semen Wash Buffer Concentrate (SEW) (DS0038) as follows:**

Number of Preps	Semen Wash Buffer Concentrate (SEW)	Molecular Biology Grade Water
20	12 ml	228 ml
50	30 ml	570 ml
250	150 ml	2850 ml

NOTE: User should carry out dilution of Semen Wash Buffer Concentrate (SEW) in a separate container, as it is not provided with this kit.

8. **Dilute Wash Solution Concentrate (WS) (DS0012) as follows:**

Number of Preps	Wash Solution Concentrate (WS)	Ethanol (96-100 %)
20	10 ml	30 ml
50	25 ml	75 ml
250	125 ml	375 ml

9. Ensure that clean & dry tubes and tips are used for the procedure.

RNase A enzyme treatment

RNase A is a type of RNase that is commonly used in research. RNase A (e.g., bovine pancreatic ribonuclease A) is one of the sturdiest enzymes in common laboratory usage. It cleaves 3'end of unpaired C and U residues.

Unit Definition for RNase A

One unit of the enzyme causes an increase in absorbance of 1.0 at 260 nm when yeast RNA is hydrolyzed at 37°C and pH 5.0. Fifty units are approximately equivalent to 1 Kunitz unit. It is completely free of DNases and proteases. The specific activity is more than 90 U/mg.

The product as supplied is stable at room temperature (15-25°C).

Centrifugation

All centrifugation steps are carried out in conventional laboratory centrifuge e.g. Beckman CS-6KR, Heraeus Varifuge 3.0R, or Sigma 6k10 with fixed angle rotor. The tubes provided with the kit are compatible with almost all laboratory centrifuges and rotors. All centrifugation steps are performed at room temperature and are given in g, the correct rpm can be calculated using the formula:

$$RPM = \sqrt{RCF/1.118} \times 10^{-5} r$$

where *RCF* = required gravitational acceleration (relative centrifugal force in units of g); *r* = radius of the rotor in cm; and *RPM* = the number of revolutions per minute required to achieve the necessary *g*-force.

Specimen Handling and Collection

Collect human/animal semen sample in a sterile container and freeze the sample at -20°C for short term storage or -80°C for long term storage. Ensure that the sample is at room temperature before beginning the protocol.

Types of Specimen

Clinical/ Animal samples: semen sample

Procedure

1. **Washing of Semen Sample**

Add 10 ml of diluted Semen Wash Buffer (SEW) (DS0038) (**Refer Step 5 of General Preparation Instructions**) to 100-200 µl of semen sample in a Corex® (or any other suitable material) centrifuge tube and vortex at full speed for 10 seconds to ensure proper mixing.

NOTE: Corex® tubes are recommended as sperm cells do not adhere strongly to the material of these tubes.

NOTE: If only Sperm DNA is required, the other cells present in the semen sample, including lymphocytes, should be separated out using appropriate gradient technique, such as

Lymphocyte Separation Medium (LSM) (Product Code LS001), to obtain only sperms in the sample which will be treated further for extraction of DNA.

2. Centrifuge at $\geq 2,500 \times g$ ($\approx 4,000$ rpm) for 10 minutes.
3. Discard the supernatant, leaving behind approximately 1 ml of pellet and Semen Wash Buffer (SEW). Vortex for 10–15 seconds to ensure proper mixing of pellet and diluted Semen Wash Buffer (SEW), and transfer the contents into a 2.0 ml capped collection tube.
4. To the above sample, add 0.5 ml of diluted Semen Wash Buffer (SEW) and vortex for 10–15 seconds to collect any sample adhering to the walls of the tube.
5. Centrifuge at $\geq 6,500 \times g$ ($\approx 10,000$ rpm) for 2 minutes. Carefully remove the supernatant, avoiding removal of the semen pellet.

6. Resuspension of pellet

Resuspend the pellet in Resuspension Solution (1X PBS) (ML116) to make the final volume 100 μ l.

7. Treatment with Sperm Lysis Buffer (SL) (DS0039)

(Prepare Sperm Lysis Buffer as indicated in General Preparation Instructions)

Add 100 μ l of prepared Sperm Lysis Buffer (SL) (DS0039) to 100 μ l of resuspended semen pellet from Step 6. Mix thoroughly by vortexing for 10-15 seconds. Incubate at 55°C for 1½ - 2 hours to ensure thorough lysis of the sample. Invert the tube intermittently for at least 8-10 times for homogenous lysis of the sample during the incubation period. If residual RNA is not a concern, continue with step 8 to treat the semen lysate further.

NOTE: More of intermittent shaking of the treated sample and less of pulse vortexing is recommended to prevent shearing of the sperm DNA.

NOTE: If reduced incubation time is desired, the user will have to optimize the incubation time at 55°C accordingly, to ensure complete homogenization of the sample.

Optional RNase A treatment

If RNA-free genomic DNA is required, add 20 μ l of RNase A Solution (DS0003) and incubate for 2 minutes at room temperature (15-25°C), then continue with step 8.

8. Lysis

Add 200 μ l of Lysis Solution (C1) (DS0010) to 200 μ l of semen lysate from step 7. Mix by vortexing thoroughly for 10-15 seconds. A homogenous mixture is essential for efficient cell lysis. Incubate at 70°C for 10 minutes.

NOTE: If any clumps are visible, the sample can be pipetted gently to obtain a homogenous mixture.

9. Prepare for Binding

Add 200 μ l of ethanol (96-100 %) to the lysate and mix thoroughly by vortexing for 5-10 seconds.

NOTE: A homogenous solution is essential.

10. Load lysate in HiElute Miniprep Spin Column (Capped) [DBCA03]

Transfer the lysate obtained from step 9 onto the column provided. Centrifuge at $\geq 6,500 \times g$ ($\approx 10,000$ rpm) for 1 minute. Discard the flow-through liquid and place the column in a same 2.0 ml uncapped collection tube.

NOTE: Use a wide bore pipette tip to reduce shearing of the DNA while transferring contents into the column.

11. First Wash

(Prepare Wash Solution as indicated in General Preparation Instructions)

Add 700 μ l of diluted Wash Solution (WS) to the column. Centrifuge at $\geq 6,500 \times g$ ($\approx 10,000$ rpm) for 1 minute. Discard the flow-through liquid and re-use the same collection tube with the column.

12. Second Wash

Add another 700 μ l of diluted Wash Solution (WS) to the column and centrifuge at 12,000-16,000 $\times g$ ($\approx 13,000$ -16,000 rpm) for 3 minutes to dry the column. Discard the flow-through liquid and centrifuge the column for another minute at the same speed if residual ethanol is observed

NOTE: (i) The column must be free of ethanol before eluting the DNA.

(ii) The collection tube can be emptied and re-used for this additional centrifugation step.

13. DNA Elution

Place the column in a new 2.0 ml uncapped collection tube. Pipette 100 μ l of the Elution Buffer (ET) (DS0040) directly onto the column without spilling to the sides. Incubate for 1 minute at room temperature (15-25°C). Centrifuge at $\geq 6,500 \times g$ ($\approx 10,000$ rpm) for 1 minute to elute the DNA. Repeat the step again with another 100 μ l of Elution Buffer (ET) for high yield of DNA.

NOTE: DNA elution can also be performed in single step by the addition of 200 μ l of Elution Buffer (ET) at a time (DNA yield would be low). Storing DNA in water may cause acid hydrolysis. To increase the elution efficiency, incubate for 5 minutes at room temperature (15-25°C) after adding the Elution Buffer (ET), then centrifuge. Elution with volume less than 200 μ l increases the final DNA concentration in the eluate significantly, but slightly reduces the overall DNA yield. Storing DNA in water may cause acid hydrolysis.

14. Transfer the eluate to a fresh capped 2ml collection tube for longer DNA storage.

Storage of the eluate with purified DNA: The eluate contains pure genomic DNA. For short-term storage (24-48 hrs) of the DNA, 2-8°C is recommended. For long-term storage, -20°C or lower temperature (-80°C) is recommended. Avoid repeated freezing and thawing of the sample which may cause denaturing of DNA. The Elution Buffer will help stabilize the DNA at these temperatures.

Alternative Short Protocol

1. Treatment with Sperm Lysis Buffer (SL)

(Prepare Sperm Lysis Buffer as indicated in General Preparation Instructions)

Add 100 μ l of prepared Sperm Lysis Buffer (SL) (DS0039) to 100 μ l of semen sample. Mix thoroughly by vortexing for 10-15 seconds. Incubate at 55°C for 1½ - 2 hours to ensure thorough lysis of the sample. Invert the tube intermittently for at least 8-10 times for homogenous lysis of the sample during the incubation period. If residual RNA is not a concern, continue with step 2 to treat the semen lysate further.

NOTE: More of intermittent shaking of the treated sample and less of pulse vortexing is recommended to prevent shearing of the sperm DNA.

NOTE: If reduced incubation time is desired, the user will have to optimize the incubation time at 55°C accordingly, to ensure complete homogenization of the sample.

Optional RNase A treatment

NOTE: RNase A treatment is generally not required while working with direct semen sample lysis method. The RNase in the seminal plasma is generally sufficient to ensure degradation of the sperm RNA.

However, if RNase treatment is still needed, add 20 µl of RNase A Solution (DS0003) to semen lysate from Step 1 and incubate for 2 minutes at room temperature (15-25°C). Continue with step 2.

2. Lysis

Add 200µl of Lysis Solution (C1) (DS0010) to 200µl of semen lysate from step 1. Mix by vortexing thoroughly for 15 seconds. A homogenous mixture is essential for efficient cell lysis. Incubate at 70°C for 10 minutes

NOTE: If any clumps are visible, the sample can be pipetted gently to obtain a homogenous mixture.

3. Continue with steps 9-13 of the regular procedure to extract sperm DNA from the sample.

Warning and Precautions

Certified for *in vitro* Diagnostic Use (IVD). Not for Medicinal Use. Read the procedure carefully before beginning the protocol. Wear protective gloves/protective clothing/eye protection/face protection. Follow good clinical laboratory practices while handling clinical samples. Standard precautions should be followed as per established guidelines. Safety guidelines may be referred in safety data sheets of the product.

Limitations

1. The yield of DNA depends upon the type and the volume of starting material used.

Performance and Evaluation

Performance of the kit is expected when the kit is used as per the protocol mentioned in the product insert within the expiry period when stored at recommended temperature.

Quality Control

Type of Sample	DNA Purity
Semen sample	1.6-1.9

References

1. Sambrook, J., *et al.* Molecular Cloning: A laboratory Manual, 2nd ed. (Cold Spring Harbor Laboratory Press, Plainview, NY, 1989)
2. Birren, B. and Lai, E. Pulsed Field Gel Electrophoresis: A practical guide (Academic Press, San Diego, CA, 1993)

Troubleshooting guide

Sr. No.	Problem	Possible Cause	Solution
1.	Spin column is clogged	Lysate/ethanol mixture is not homogenous	Vortex the tubes for atleast 5-10 seconds in order to obtain a homogenous solution before applying it to the column. If minimally sheared DNA is required for downstream applications like PCR, mix with gentle pipetting or inversion until homogenous lysate is obtained, instead of vortexing.
		DNA elution is improper	Ensure that the DNA elution is in 200µl of Elution Buffer. To improve the DNA yield incubate for 5 minutes at room temperature after Elution Buffer is added to the column.
		Ethanol was omitted during binding	Ensure that ethanol is added in step 9 before adding the sample to the spin column in step 10.
		Eluate contains residual ethanol from the wash	Remove ethanol from the second wash completely before eluting the DNA. Spin for an additional 2 minute to dry the membrane completely. In order to avoid the interference of ethanol, always use a fresh tube for elution.
		Use of water instead of Elution Buffer for elution of DNA	Elution Buffer is recommended for optimal yields and storage of the genomic DNA. If water is used instead of the Elution Buffer the pH should be at least 7.0, to avoid acidic conditions which may cause acid hydrolysis of DNA when stored for long periods of time. NOTE: Only DNase/RNase and Protease free water should be used for eluting DNA.
2.	Purity of the DNA is lower than expected; (A_{260}/A_{280} ratio is low)	Eluate was diluted in water for absorbance measurement.	Use either the Elution Buffer provided, or 10 mM Tris-HCl, pH 8.5.
		Background reading is high due to silica fines	Spin the DNA sample at maximum speed for 1 minute, the supernatant can be used to repeat the absorbance readings
3.	Purity of the DNA is higher than expected; (A_{260}/A_{280} ratio is too high)	RNA contamination	RNase A treatment can be included in future isolations or the final product can be treated with RNase A Solution and repurified.

4.	Shearing of genomic DNA	Improper handling of genomic DNA	All pipetting steps should be carried as gently as possible. Wide bore pipette tips are recommended to eliminate shearing of the DNA to a large extent. If the isolated DNA is to be used for PCR, mix with gentle pipetting or invert until homogenous solution is obtained to reduce shearing of DNA considerably.
5.	Downstream applications are inhibited	Traces of ethanol present in the final genomic DNA preparation	After the washing steps the eluate should not come in contact with the column. Spin the column for 1 minute at maximum speed (12,000-16,000 x g) if necessary, after emptying the collection tube.
		Salt is carried over in the final genomic DNA preparation	The spin column should be transferred to a new 2.0 ml collection tube before adding the elution buffer.

Safety Information



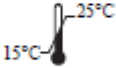


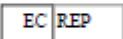

The HiPurA® Sperm Genomic DNA Purification Kit is for laboratory use only, not for drug, household or other uses. The Lysis Solution (C1) contains chaotropic salts, which are irritant. Take appropriate laboratory safety measures and wear gloves when handling. Not compatible with disinfecting agents containing bleach.

Disposal

User must ensure safe disposal by autoclaving and/or incineration of used or unusable preparations of this product. Follow established laboratory procedures in disposing of infectious materials and material that comes into contact with clinical sample must be decontaminated and disposed off in accordance with current laboratory techniques.

Technical Assistance

At HiMedia, we pride ourselves on the quality and availability of our technical support. For any kind of technical assistance, send an email to mb@himedialabs.com.

	In vitro diagnostic medical device
	CE Marking
	Storage temperature
	Do not use if package is damaged
	HiMedia Laboratories Private Limited, Reg. Off: Plot No. C-40, Road No. 21Y, MIDC, Wagle Industrial Area, Thane, (West) 400604, Maharashtra, INDIA. Web: www.himedialabs.com
	CE Partner 4U ,Esdoornlaan 13, 3951 DB Maarn The Netherlands, www.cepartner4u.eu
	02/2025

PIMB522_0/0222

MB522-09

Disclaimer :

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal or therapeutic use but for laboratory, diagnostic, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.

HiMedia Laboratories Pvt. Ltd. Reg.office : Plot No. C-40, Road No. 21Y, MIDC, Wagle Industrial Area, Thane, (West) 400604, Maharashtra, INDIA.
Customer Care No.: 00-91-22-6116 9797 Tel: 00-91-22-6147 1919, 6903 4800 Email: techhelp@himedialabs.com Website: www.himedialabs.com