

MBT075

Hi-PCR Kit

Product Name	Product Code	Kit Packing
Hi-PCR Kit	MBT075-10R	10 reactions
Hi-PCR Kit	MBT075-25R	25 reactions

Description:

Hi-PCR Kit is provided at affordable price and is designed to meet various needs while providing convenience in many ways. Hi-PCR kit contains all necessary components for PCR. The user needs to add the desired template and primers of interest. This kit is beneficial when sharing the reagents with other users.

Hi-PCR kit is provided with:

Components	Product Code	Recommended volume to be added for 50 µl reaction	Final concentration in the reaction mixture	Reagents provided for 10R (reactions)	Reagents provided for 25R (reactions)
10X Assay Buffer	DS0278	5 µl	1X	75 µl	150 µl
10 mM dNTP mix	MBT078	5 µl	1 mM	75 µl	150 µl
50 mM MgCl ₂	DS0118	2.5 µl	2.5 mM	75 µl	150 µl
Taq Polymerase (1 U/µl)	MBT060D	2.5 µl	1 Units	50 µl	100 µl
Molecular Biology Grade Water for PCR	ML065	Upto 50 µl	-	1 ml	2 ml
6X Loading Dye	ML015	-	-	100 µl	200 µl

Recommended PCR program:

1. Initial denaturation : 94°C for 5-10 minutes
2. Cycling Parameters (No. of cycles: 27-35)
 - Denaturation : 94°C for 45 seconds
 - Annealing : 50-60°C for 30 seconds
 - Extension : 72°C for 30 seconds
3. Final Extension : 72°C for 5-10 minutes

Storage conditions: Store between -10°C to -20°C. When stored under recommended conditions, the kit components are stable for 1 year.


Quality control:

Detected free of RNases, endonuclease and exonuclease activities.

Please refer disclaimer Overleaf.

Technical Assistance

At HiMedia, we pride ourselves on the quality and availability of our technical support. For any kind of technical assistance, mail at mb@himedialabs.com.

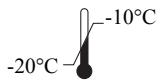

Storage temperature



Do not use if package is damaged



HiMedia Laboratories Private Limited,
Reg. Off: Plot No. C-40, Road No. 21Y,
MIDC, Wagle Industrial Area, Thane,
(West) 400604, Maharashtra, INDIA.
www.himedialabs.com



01/2025

PIMBT075_0/0122

MBT075-06

Disclaimer :

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal or therapeutic use but for laboratory, diagnostic, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.

HiMedia Laboratories Pvt. Ltd. Reg.office : Plot No. C-40, Road No. 21Y, MIDC, Wagle Industrial Area, Thane, (West) 400604, Maharashtra, INDIA.
Customer Care No.: 00-91-22-6116 9797 Tel: 00-91-22-6147 1919, 6903 4800 Email: techhelp@himedialabs.com Website: www.himedialabs.com