



## Carbol Fuchsin (ZN,Strong)

**S005**

### Intended use

Carbol Fuchsin (Zn,Strong) is used as staining solution for acid fast staining.

### Composition\*\*

#### Ingredients

Basic Fuchsin	0.30 gm
Ethyl alcohol, 95%	10.0 ml
Phenol	5.0 gm

### Directions

1. Prepare a smear on a clear, dry glass slide.
2. Allow it to air dry and fix with gentle heat.
3. Flood the smear with Carbol Fuchsin stain (S005). Heat to steaming for 5 minutes with a low flame; do not boil the stain and do not permit drying of the smear.
4. Allow it to stand for 5 minutes without further heating.
5. Wash in the running tap water.
6. Decolourize with Acid Fast Decolourizer (S033) for 2 minutes or until no more stain comes off in the washings. (If washing is not thorough, may get false positive results).
7. Wash with the tap water.
8. Counterstain for 30 seconds with Methylene Blue (S022).
9. Wash with tap water, dry in air, then examine under an oil immersion objective.

### Principle And Interpretation

Mycobacteria (AFB/Acid Fast Bacteria) are difficult to stain due to high lipid and wax content in their cell walls. This differential staining technique is useful for identification of the tubercle bacillus, other Mycobacteria and Nocardia, which depends on the chemical composition of the bacterial cell wall. Because of the difficulty in staining these organisms with ordinary dyes, basic dyes in the presence of controlled amounts of acid are used. Generally, heat must be applied during the staining procedure, or wetting agents must be used, to aid dye penetration. Organisms exhibiting the property of acid-fastness, once stained, are not easily decolourized by alcohol.

### Type of specimen

Clinical specimen: sputum

### Specimen Collection and Handling

All testing for acid-fast bacilli is sent to the reference laboratory in an effort to meet the 24 hr TAT time for smear results. Use sterile, leak proof disposable plastic containers for collection. Do not use wax containers as these can cause false positive smear results. Do not use any fixative or preservatives. Swabs are not recommended as a collection device for the isolation of mycobacteria. They are acceptable only if the specimen can not be obtained by

any other means. A negative result from a swab specimen is not reliable. In general, the number of acid fast bacilli in a specimen is small. Early morning specimens are the specimens of choice for urine and sputum because the mycobacteria have had a chance to pool and concentrate, and so increase the chances of recovery. Always collect and submit the maximum volume possible of specimens normally considered sterile. Do not submit 24-hour collections, as they are likely to be diluted and contaminated. Collect specimens before antimicrobial therapy is started. Even a few days of therapy may kill or inhibit sufficient numbers of mycobacteria to prevent recovery on culture and so leave confirmation of disease in doubt. If a specimen is submitted after therapy has been initiated, note on the request. Avoid contamination of the specimen with tap water, as environmental mycobacteria exist and their recovery by smear or culture can cause confusion for the patient diagnosis.

## Warning and Precautions

In Vitro diagnostic use only. Read the label before opening the container. Wear protective gloves/protective clothing/ eye protection/face protection. Follow good microbiological lab practices while handling specimens and culture. Standard precautions as per established guidelines should be followed while handling clinical specimens. Safety guidelines may be referred in individual safety data sheets.

## Limitations

1. Excessive washing following the carbol fuchsin may cause a heightens decolourization effect.

## Performance and Evaluation

Performance of the medium is expected when used as per the direction on the label within the expiry period when stored at recommended temperature

## Quality Control

- **Appearance** : Dark pinkish red solution, clear solution with sheen.
- **Clarity** : Clear without any particles.(Note: On storage of stain, Precipitate may develop. This will not affect the performance criteria of the stain.)
- **Microscopic Examination** : Acid fast staining (Ziehl Neelsen`s method) is carried out where Carbol Fuchsin is used as one of the stains and staining characteristic of orga nisms is observed under microscope by using oil immersion lens.
- **Results** : Acid fast organisms: Bright red  
All other organisms and background material: Blue

## Storage and Shelf Life

Store between 10-30°C in tightly closed container and away from bright light. Use before expiry date on label. On opening, product should be properly stored in dry ventilated area protected from extremes of temperature and sources of ignition. Seal the container tightly after use.

## Disposal

User must ensure safe disposal by autoclaving and/or incineration of used or unusable preparations of this product. Follow established laboratory procedures in disposing of infectious materials and material that comes into contact with clinical sample must be decontaminated and disposed of in accordance with current laboratory techniques.

## Reference

1. Downes F. P. and Ito K. (Ed.), 2001, Compendium of Methods for the Microbiological Examination of Foods,

4th ed., APHA, Washington, D.C.

2. Rice E.W., Baird, R.B., Eaton A. D., Clesceri L. S. (Eds.), 2012, Standard Methods for the Examination of Water and Wastewater, 22nd ed., APHA, Washington, D.C.

3. Wehr H. M. and Frank J. H., 2004, Standard Methods for the Microbiological Examination of Dairy Products, 17th Ed., APHA Inc., Washington, D.C.

4. Isenberg, H.D. Clinical Microbiology Procedures Handbook. 2nd Edition.

5. Jorgensen, J.H., Pfaller, M.A., Carroll, K.C., Funke, G., Landry, M.L., Richter, S.S and Warnock, D.W. (2015) Manual of

Clinical Microbiology, 11th Edition. Vol. 1.

6. Shanhoodtzer, C.J., P. Schaper, and L.R. Peterson. 1982. Concentrated Gram stain smear prepared with a cytospin centrifuge. J.clin. Microbiol.16:1052-1056

7. Thorpe, J.E., R.P. Banghman, P.T. Frame, T.A. Wessler, and J.L. Staneck. 1987. Bronchoalveolar lavage for diagnosing acute bacterial pneumonia. J.Infect.Dis.155:855-861

8. Clinical Microbiology Procedures Handbook, 1992, Isenberg, American Society for Microbiology. LabCorp Specimen Collection Instructions



Storage temperature



Do not use if package is damaged



In vitro diagnostic medical device



CE Marking



HiMedia Laboratories Pvt Limited  
C-40, 21/Y, MIDC, Wagle Ind Area,  
Thane(W)-400604, Maharashtra, India



CEpartner4U, ESDOORNLAAN 13, 3951  
DB MAARN, The Netherlands,  
[www.cepartner4u.eu](http://www.cepartner4u.eu)

Revision 02/2022

#### Disclaimer :

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal or therapeutic use but for laboratory, diagnostic, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.